

RiboCluster Profiler™

RIP-Certified Antibody

Anti-PCBP2 pAb

Code No.	Quantity	Concentration	Form
RN025P	200 µL	1 mg/mL	Affinity Purified

BACKGROUND: Poly(rC)-binding proteins (PCBPs) are cellular RNA-binding proteins, characterized by their triple KH structure and by their poly(C)-binding specificity. The 3 KH domains are not only involved in nucleic acid binding but also mediate protein-protein interactions within the RNP complexes. It binds specifically to polypyrimidine tracts within the 3' UTR of target mRNAs through KH domains. PCBPs are expressed as 4 isoforms in cells. PCBP1 and PCBP2 are predominantly expressed in the nucleus and can shuttle from the nucleus to the cytoplasm, while PCBP3 and PCBP4 are expressed to a less extent in the cytoplasm. The most characterized gene regulated by PCBP1 and PCBP2 is the α -globin gene, and the accumulated α -globin mRNA is known to account for approximately 95% of total cellular mRNA during terminal erythroid differentiation. Specific binding of PCBPs to 3' UTR of mRNA is associated with the stabilization of α -globin mRNA. Microarray analysis of human mRNAs present in a specific subset of PCBP2-RNP complexes also showed a potential involvement of PCBP2 in transcription factor expression via association with their mRNA; further, the enrichment of PCBP2 mRNA in PCBP2-RNP complexes suggested that PCBP2 expression may be subject to posttranscriptional autoregulation.

RIP-CERTIFIED ANTIBODY:

Posttranscriptional regulation of gene expression is a ribonucleoprotein-driven process, which involves RNA binding proteins (RBPs) and non-coding RNAs that affect splicing, nuclear export, subcellular localization, mRNA decay and translation. The RNP Immunoprecipitation-Chip (RIP-Chip), RIP-Seq and RIP-RT-PCR allow the identification of multiple RNA targets of RBPs globally and within the context of a cell extract. Antibodies specific to the RNA binding protein of interest are used to co-immunoprecipitate the RNA binding protein and the associated subset of mRNAs. The mRNA content is interrogated using standard microarray or sequencing technology. RIP-Certified Antibody is validated for use in RNP Immunoprecipitation (RIP) in conjunction with the RIP-Assay Kit distributed from MBL. Its ability to immunoprecipitate mRNAs and RBPs complex was confirmed by quantitative and qualitative analysis on NanoDrop, Bioanalyzer and RT-PCR or microarray.

SOURCE: This antibody was purified from rabbit serum by affinity column chromatography. The rabbit was immunized with KLH conjugated synthetic peptide, corresponding to internal region of human PCBP2.

FORMULATION: 200 µL volume of PBS containing 50% glycerol, pH 7.2. No preservative is contained.

STORAGE: This antibody solution is stable for one year from the date of purchase when stored at -20°C.

REACTIVITY: This antibody reacts with human PCBP2 on Western blotting, Immunoprecipitation and RNP Immunoprecipitation.

INTENDED USE:

For Research Use Only. Not for use in diagnostic procedures.

APPLICATIONS:

RNP Immunoprecipitation: 15 µg/500 µL of cell extract from 1.5×10^7 cells

Western blotting: 1 µg/mL for chemiluminescence detection system

Immunoprecipitation: 5 µg/500 µL of cell extract from 5×10^6 cells

Immunohistochemistry: Not tested

Immunocytochemistry: Not tested*

Flow cytometry: Not tested

Crosslinking-immunoprecipitation (CLIP): Not tested*

*It is reported that this antibody can be used in Immunocytochemistry^{1), 3)}, enhanced CLIP^{2), 3)} and Immunodepletion⁵⁾.

Detailed procedures are provided in the following **PROTOCOLS**.



SPECIES CROSS REACTIVITY:

Species	Human	Mouse	Rat	Hamster
Cells	293T, HeLa, K562, Jurkat, MCF7	NIH/3T3, WR19L	Rat1	Not tested
Reactivity on WB	+	+	+	

REFERENCES:

- 1) Markmiller, S., *et al.*, *Cell* **172**, 590-604.e13 (2018) [IC]
- 2) Van Nostrand, E. L., *et al.*, *Nat. Methods* **13**, 508-514 (2016) [CLIP]
- 3) Sundararaman, B., *et al.*, *Mol. Cell* **61**, 903-913 (2016) [IC, CLIP]
- 4) Neff, A. T., *et al.*, *Genome Res.* **22**, 1457-1467 (2012) [WB]
- 5) Palusa, S., *et al.*, *PLoS One* **7**, e33561 (2012) [RIP]
- 6) Waggoner, S. A., and Liebhaber, S. A., *Mol. Cell. Biol.* **23**, 7055-7067 (2003)

- 7) Chkheidze, A. N., and Liebhaber, S. A., *Mol. Cell. Biol.* **23**, 8405-8415 (2003)
8) Makeyev, A. V., and Liebhaber, S. A., *RNA* **8**, 265-278 (2002)

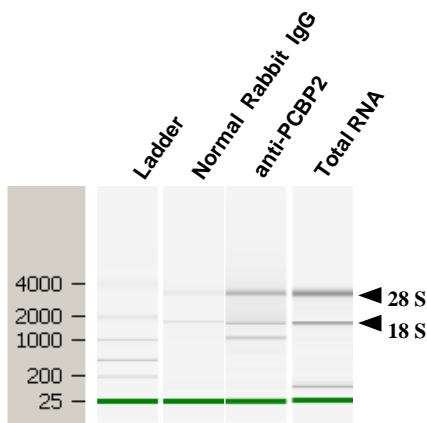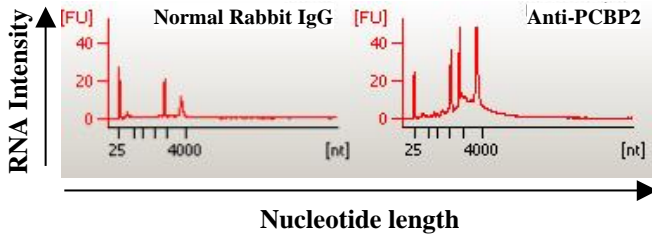

Analysis of isolated RNA with Bioanalyzer.

Average of the RNA Quantity (n=2)	
Antibody	RNA (ng)
Normal Rabbit IgG	89.0
anti-PCBP2	369.0
Total RNA	230260.0

PROTOCOLS:

RNP Immunoprecipitation

Some buffers and reagents are included in the RIP-Assay Kit (code. RN1001). Please also refer to the protocol packaged in the RIP-Assay Kit.

[Material Preparation]

1. Lysis Buffer (+)
Before using the Lysis Buffer, protease inhibitors, RNase inhibitors, and DTT are added to the Lysis Buffer at the appropriate concentration.
2. Wash Buffer (+)
Before using the Wash Buffer, DTT is added to the Wash Buffer at the appropriate concentration.


Protocol

- 1) Wash 1.5×10^7 cells 4 times with PBS and resuspend them with 500 μ L of ice-cold Lysis Buffer (+) containing appropriate protease inhibitors, RNase inhibitors, and DTT. Vortex thoroughly, then incubate it on ice for 10 minutes.
- 2) Centrifuge the tube at 12,000 x g for 5 minutes at 4°C and transfer the supernatant to another tube.
- 3) Add 25 μ L of 50% protein A agarose beads slurry

resuspended in Lysis Buffer (+) into the supernatant. Incubate it at 4°C with rotating for 1 hour.

- 4) Centrifuge the tube at 2,000 x g for 1 minute at 4°C and transfer the supernatant to another fresh tube (precleared sample).
- 5) Mix 25 μ L of 50% protein A agarose beads slurry resuspended in nuclease-free PBS with Normal Rabbit IgG (RIP-Assay Kit) or Anti-PCBP2 pAb (RN025P) at the amount as suggested in the **APPLICATIONS**, and then add 1 mL of Wash buffer (+) into each tube. Incubate with gentle agitation for 1 hour at 4°C.
- 6) Wash the beads once with ice-cold Lysis Buffer (+) (centrifuge the tube at 2,000 x g for 1 minute). Carefully discard the supernatant using a pipettor without disturbing the beads.
- 7) Add 500 μ L of cell lysate (precleared sample of step 4), then incubate with gentle agitation for 3 hours at 4°C.
- 8) Wash the bead pellet 4 times with Wash Buffer (+) (centrifuge the tube at 2,000 x g for 1 minute).
- 9) Add 400 μ L of Master mix solution (Solution I: Solution II = 10 μ L: 390 μ L). Vortex thoroughly, then spin-down.
- 10) Add 250 μ L of Solution III. Vortex thoroughly.
- 11) Centrifuge the tube at 2,000 x g for 2 minutes.
- 12) Transfer the supernatant to the fresh tube containing 2 μ L of Solution IV.
- 13) Add 600 μ L of ice-cold 2-propanol and place at -20°C for 20 minutes. Centrifuge the tube at 12,000 x g for 10 minutes.
- 14) Wash the pellet 2 times with 0.5 mL of ice-cold 70% Ethanol and dry up the pellet for 5-15 minutes.
- 15) Dissolve the pellets in nuclease-free water.
- 16) RNA was quantified with NanoDrop (Thermo Fisher Scientific Inc.) and the RNA quality was analyzed with Bioanalyzer (Agilent Technologies, Inc.).

(Positive control for RNP Immunoprecipitation; K562)

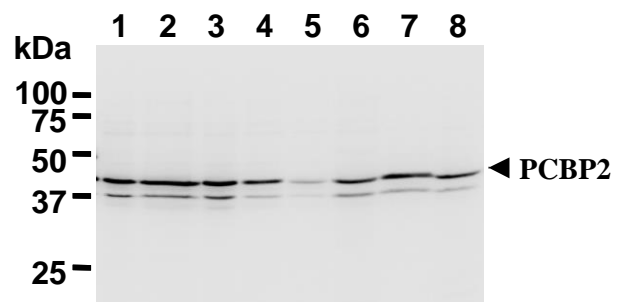


Western blot analysis of PCBP2 expression in 293T (1), HeLa (2), K562 (3), Jurkat (4), MCF7 (5), NIH/3T3 (6) WR19L (7) and Rat1 (8) using RN025P.

SDS-PAGE & Western Blotting

- 1) Wash 1×10^7 cells 3 times with PBS and suspend them in 1 mL of Laemmli's sample buffer.
- 2) Boil the samples for 2 minutes and centrifuge. Load 10 μ L of sample per lane on a 1-mm-thick SDS-polyacrylamide gel and carry out electrophoresis.

- 3) Blot the protein to a polyvinylidene difluoride (PVDF) membrane at 1 mA/cm² for 1 hour in a semi-dry transfer system (Transfer Buffer: 25 mM Tris, 190 mM glycine, 20% MeOH). See the manufacturer's manual for precise transfer procedure.
- 4) To reduce nonspecific binding, soak the membrane in 10% skimmed milk (in PBS, pH 7.2) for 1 hour at room temperature, or overnight at 4°C.
- 5) Incubate the membrane with primary antibody diluted with 1% skimmed milk (in PBS, pH 7.2) as suggested in the **APPLICATIONS** for 1 hour at room temperature. (The concentration of antibody will depend on the conditions.)
- 6) Wash the membrane with PBS-T [0.05% Tween-20 in PBS] (5 minutes x 3 times).
- 7) Incubate the membrane with 1:10,000 of Anti-IgG (Rabbit) pAb-HRP (MBL; code no. 458) diluted with 1% skimmed milk (in PBS, pH 7.2) for 1 hour at room temperature.
- 8) Wash the membrane with PBS-T (5 minutes x 3 times).
- 9) Wipe excess buffer off the membrane, and incubate membrane with an appropriate chemiluminescence reagent for 1 minute.
- 10) Remove extra reagent from the membrane by dabbing with a paper towel, and seal it in plastic wrap.
- 11) The detection was performed with LAS-4000 (FUJIFILM).

(Positive controls for Western blotting; 293T, HeLa, K562, Jurkat, MCF7, NIH3T3, WR19L and Rat1)

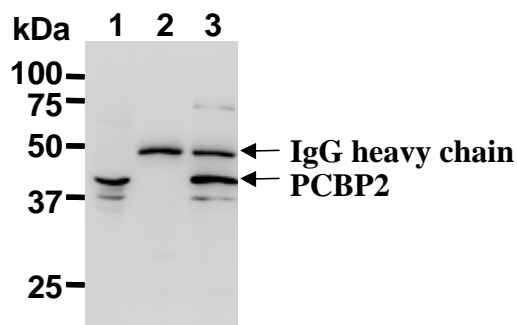

Immunoprecipitation of PCBP2 from K562 with normal rabbit IgG (2) or RN025P (3). After immunoprecipitated with the antibody, immunocomplex was resolved on SDS-PAGE and immunoblotted with RN025P. Lane 1 is the input sample.

Immunoprecipitation

- 1) Wash cells (approximately 1 x 10⁷ cells) 2 times with PBS and resuspend them with 1 mL of ice-cold Lysis buffer (RIP-Assay Kit) containing protease inhibitors and DTT at appropriate concentrations. Vortex for 10 seconds. Leave on ice for 10 minutes.
- 2) Centrifuge the tube at 12,000 x g for 5 minutes at 4°C and transfer the supernatant to another tube.
- 3) Add 20 µL of 50% protein A agarose beads slurry resuspended in Lysis Buffer into the supernatant. Incubate it at 4°C with rotating for 1 hour.
- 4) Centrifuge the tube at 2,000 x g for 1 minute at 4°C and

transfer the supernatant to another tube (precleared sample).

- 5) Mix 20 µL of 50% protein A agarose beads slurry resuspended in PBS with Normal Rabbit IgG (RIP-Assay Kit) or Anti-PCBP2 pAb (RN025P) at the amount as suggested in the **APPLICATIONS**, and then add 1 mL of Wash buffer into each tube. Incubate with gentle agitation for 1 hour at 4°C.
- 6) Wash the beads once with ice-cold Lysis Buffer (centrifuge the tube at 2,000 x g for 1 minute). Carefully discard the supernatant using a pipettor without disturbing the beads.
- 7) Add 500 µL of cell lysate (precleared sample of step 4), then incubate with gentle agitation for 3 hours at 4°C.
- 8) Wash the beads 4 times with Wash Buffer (centrifuge the tube at 2,000 x g for 1 minute).
- 9) Resuspend the beads in 20 µL of Laemmli's sample buffer, boil for 3-5 minutes, and centrifuge for 5 minutes. Use 20 µL/lane for the SDS-PAGE analysis. (See **SDS-PAGE & Western blotting.**)

(Positive control for Immunoprecipitation; K562)

RELATED PRODUCTS:

RIP-Certified Antibody

RN001M	Anti-IGF2BP1 (IMP1) mAb (6H6)
RN004M	Anti-Ribosomal P0/P1/P2 mAb (9D5)
RN005M	Anti-EIF2C2 (AGO2) mAb (2A8)
RN006M	Anti-EIF4E mAb (C107-3-5)
RN007M	Anti-ELAVL1 (HuR) mAb (C67-1)
RN009M	Anti-PABPC1 mAb (10E10)
RN001P	Anti-EIF4E pAb
RN002P	Anti-EIF4G1 (Human) pAb
RN003P	Anti-EIF4G2 pAb
RN004P	Anti-ELAVL1 (HuR) pAb
RN005P	Anti-ELAVL2 (HuB) (Human) pAb
RN006P	Anti-ELAVL3 (HuC) pAb
RN007P	Anti-IGF2BP1 (IMP1) pAb
RN008P	Anti-IGF2BP2 (IMP2) pAb
RN009P	Anti-IGF2BP3 (IMP3) pAb
RN010P	Anti-MSI1 (Musashi1) pAb
RN011P	Anti-PTBP1 (Human) pAb
RN012P	Anti-STAU1 (Human) pAb
RN013P	Anti-STAU2 (Human) pAb
RN014P	Anti-TIA1 pAb
RN015P	Anti-YBX1 pAb
RN016P	Anti-FMR1 pAb
RN017P	Anti-FXR1 pAb
RN018P	Anti-FXR2 pAb
RN019P	Anti-HNRNPK pAb
RN020P	Anti-ILF3 (Human) pAb
RN021P	Anti-KHDRBS1 pAb
RN022P	Anti-PABPC4 pAb
RN024P	Anti-PCBP1 pAb
RN026P	Anti-PUM1 pAb
RN027P	Anti-PUM2 pAb
RN028P	Anti-EIF2C1 (AGO1) pAb
RN032P	Anti-CIRBP pAb
RN033P	Anti-TNRC6A (GW182) (Human) pAb
RN037P	Anti-AUH pAb

RN038P Anti-CPEB1 pAb
RN041P Anti-KHDRBS2 (SLM1) pAb
RN045P Anti-SLBP pAb

RBP Antibody

RN002MW Anti-CUGBP1 mAb (3B1)
RN008MW Anti-ELAVL1 (HuR) mAb (C54-6)
RN010MW Anti-PIWIL1 (MIWI) mAb (2D9)
RN012MW Anti-Coil (Coilin) mAb (#4)
RN013MW Anti-Nono (P54NRB) mAb (C5)
RN014MW Anti-SFPQ (PSF) mAb (C23)
RN015MW Anti-PSPC1 (PSP1) mAb (1L4)
RN023PW Anti-PABPN1 pAb
RN028PW Anti-EIF2C1 (AGO1) pAb
RN029PW Anti-EIF2C2 (AGO2) pAb
RN030PW Anti-DICER1 (Human) pAb
RN031PW Anti-ZFP36 (Human) pAb
RN034PW Anti-CUGBP1 pAb
RN035PW Anti-CUGBP2 pAb
RN036PW Anti-ACO1 (IRP1) pAb
RN039PW Anti-CPEB2 pAb
RN040PW Anti-CPEB4 pAb
RN042PW Anti-MBNL1 pAb
RN043PW Anti-NOVA1 (Human) pAb
RN044PW Anti-NOVA2 (Human) pAb
RN046PW Anti-SYNERIP (HNRNPQ) pAb
RN047PW Anti-PTBP2 pAb
RN048PW Anti-G3BP1 pAb
RN049PW Anti-G3BP2 pAb
RN050PW Anti-GRSF1 pAb
RN051PW Anti-HDLBP (Vigilin) pAb
RN052PW Anti-HNRNPC pAb
RN053PW Anti-PAIP1 pAb
RN054PW Anti-PCBP3 pAb
RN055PW Anti-AIMP1 (SCYE1) pAb
RN056PW Anti-SERBP1 pAb
RN057PW Anti-TARBP1 pAb
RN058PW Anti-TARBP2 pAb
RN059PW Anti-TIAL1 pAb
RN060PW Anti-HNRNPD (AUF1) pAb
RN061PW Anti-HNRNPA0 pAb
RN062PW Anti-DGCR8 pAb
RN063PW Anti-DHX9 pAb
RN064PW Anti-FUSIP1 (SRSF10) pAb
RN065PW Anti-KHSRP pAb
RN066PW Anti-KIAA0020 pAb
RN067PW Anti-PPP1R10 pAb
RN068PW Anti-PPP1R8 pAb
RN069PW Anti-RBM14 pAb
RN070PW Anti-RPS10 pAb
RN071PW Anti-RPS19 pAb
RN072PW Anti-RPS6 pAb
RN073PW Anti-RPS9 pAb
RN074PW Anti-SSB (La) pAb
RN075PW Anti-PPARGC1B pAb
RN076PW Anti-PPRC1 pAb
RN077PW Anti-SMN1 pAb
RN078PW Anti-SMNDC1 pAb
RN079PW Anti-SRSF7 (9G8) pAb
RN080PW Anti-SRSF3 (SRp20) pAb
RN081PW Anti-SRSF9 (SRp30c) pAb

RN082PW Anti-SRSF5 (SRP40) pAb
RN083PW Anti-AQ (IBP160) pAb
RN084PW Anti-SRRM1 (SRM160) pAb
RN085PW Anti-U2AF1 pAb
RN086PW Anti-U2AF2 pAb
RN087PW Anti-ALYREF (THOC4) pAb
RN088PW Anti-NXF1 (TAP) pAb
RN089PW Anti-MAGOH pAb
RN090PW Anti-DDX21 pAb
RN091PW Anti-DDX23 pAb
RN092PW Anti-NONO (P54NRB) pAb
RN093PW Anti-PRPF4 pAb
RN094PW Anti-PRPF8 pAb
RN095PW Anti-SNRNP200 pAb
RN096PW Anti-SNRNP40 pAb
RN097PW Anti-SNRNP70 pAb
RN098PW Anti-EDC4 pAb
RN099PW Anti-EIF4A1 pAb
RN100PW Anti-EXOSC5 (RRP46) (Human) pAb
RN101PW Anti-FBL (Fibrillarin) pAb
RN102PW Anti-GEMIN2 (Human) pAb
RN103PW Anti-NCBP1 (CBP80) pAb
RN104PW Anti-PAN2 (USP52) (Human) pAb
RN105PW Anti-PARN pAb
RN106PW Anti-SFPQ (PSF) pAb
RN107PW Anti-TARDBP (TDP-43) pAb
RN108PW Anti-UPF1 pAb
RN109PW Anti-XRN1 (Human) pAb
RN110PW Anti-CNOT7 (CAF1) pAb
RN111PW Anti-ETF1 (eRF1) pAb
RN112PW Anti-DCP1B (Human) pAb
RN113PW Anti-DHX36 (RHAU) pAb
RN114PW Anti-HNRNPA1 pAb
RN115PW Anti-LIN28B (Human) pAb
RN116PW Anti-DDX39B (UAP56) pAb
RN117PW Anti-CCAR2 (DBC1) pAb
RN118PW Anti-UPF3B pAb
RN119PW Anti-GSPT2 (eRF3b) (Human) pAb
RN120PW Anti-RBM8A (Y14) pAb
RN121PW Anti-FTO (Human) pAb
RN122PW Anti-ALKBH5 pAb
RN123PW Anti-YTHDF2 pAb
RN124PW Anti-RNMT (Human) pAb
RN125PW Anti-HENMT1 pAb
RN126PW Anti-AHCY (SAHH) pAb
RN127PW Anti-NSUN2 (Human) pAb
RN128PW Anti-TRMT6 (Human) pAb
RN129PW Anti-DDX6 (RCK/p54) pAb
RN130PW Anti-TRMT61A (Human) pAb

Kit, others (reagents, equipments, etc.)

RN1001 RIP-Assay Kit
RN1005 RIP-Assay Kit for *microRNA*
RN1007 BRIC Kit
RN1011 RiboTrap Kit
MJS002V2 Protein G-Magnetic Beads
3190 Magnetic Rack

For the latest information of RiboCluster Profiler™, please visit our website at <http://ruo.mbl.co.jp/je/rip-assay/>